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510006M01

Introduction

Thank you for purchasing hanger fitting for the FMA series.

This is for under-weighing used for FMA series.


Read this instruction carefully.

 CAUTION	Due to the heavy weight of this product, lifting and carrying during assembly and installation must be carried out by two or more persons, according to the operation manual of the FMA series. Follow this manual for the installation of the hanger fitting option.
--------------------------------------------------------------------------------------------------	--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------

Note

- When a hanger fitting is installed, do not place the scale on the flat surface. The fitting may be broken, or the scale is damaged.
- When a hanger fitting is installed, the weighing unit does not meet IP65.
Do not expose the scale to liquid nor dust. Liquid or dust may infiltrate into the weighing unit and may cause to malfunction of the scale.
- This option cannot be used for legal for trade purposes.


1. Bundled Items

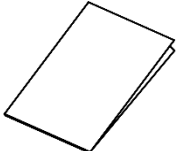1 FMA hanger fitting: 1 	2 Spacer: 1 	3 Operation manual: 1 
----------------------------------------------------------------------------------------------------------------	----------------------------------------------------------------------------------------------------	----------------------------------------------------------------------------------------------------------------


2. Installation

① Up-side down the FMA weighing unit


For scale with pole (pole type), remove the pole before up-side down the unit.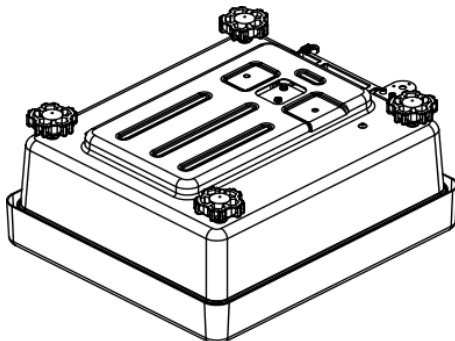

② Remove the under-weighing cover

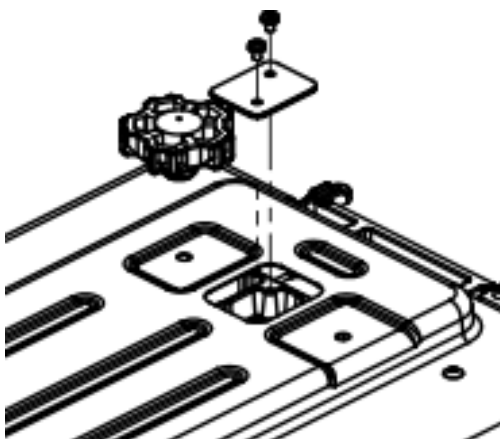③ Put a spacer and install the hanger fitting by screwing.

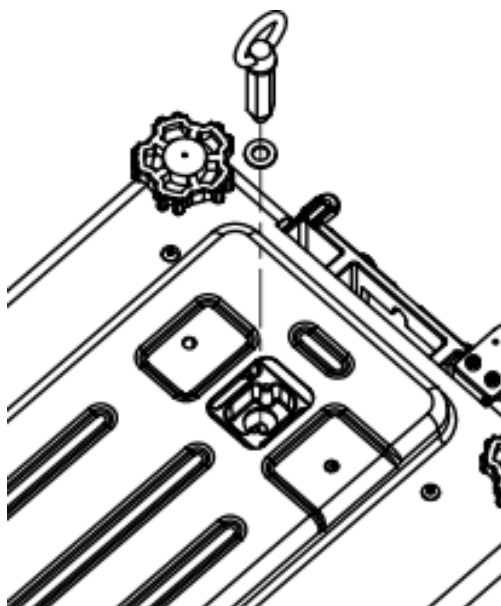3. Adjusting under-weighing pan

- ① Hang an under-weighing pan on the hanger fitting.
- ② Turn on power of the scale.
- ③ When <Net> is displayed (For non-verified scale only)
 - This indicates that the weighing range is reduced and cannot weigh up to the full capacity. Reduce the (total) weight of weighing pan(s) by using lighter under-weighing pan or by removing weighing pan of the scale.
- ④ When <ERROR 706> occurs
 - The under-weighing pan is too heavy or too light. Use lighter or heavier under-weighing pan, or replace the weighing pan of the scale.

4. Span adjustment

When usually using with the under-weighing pan, execute span adjustment using the under-weighing pan. But, when <Net> is displayed in the above “3. Adjusting under-weighing pan”, remark the following A & B when span adjustment/span test (see operation manual for the scale) is performed.

A. Semi-automatic span adjustment with internal weight, Span test with internal/external weight

If <ERR 719> appears, span adjustment/span test is not possible because the under-weighing pan is too heavy. Do such adjustment or test by removing the under-weighing pan. Alternatively, execute span adjustment by external weight.

B. When span adjustment in performed with external weight

When <Net> is displayed, weighing range is reduced and external weight of which weight is equal to the maximum capacity of the scale cannot be used. In this case, use external weight that do not cause <OVER ERROR>, select <VAR> at <SELECT MIN> and enter the mass of the external weight to execute span adjustment with external weight.